

# Dip Sampler, plunging siphon



Dip sampler for simple removal of liquid samples. Good view due to transparent tubes. Upper end with reflux valve, lower end with ball valve.

- PP, transparent
- Dip sampler, item no. 5315-0200, steplessly extendable from 100 cm to 200 cm
- All-layer and point sampling possible



# Use

## All-layer sample

- Insert the dip sampler slowly into the liquid with the valve open.
- 2. Close valve.
- 3. Pull the dip sampler back and take out the sample.

### Point sampling/bottom sample

- Insert the dip sampler to the targeted point/to the ground with closed valve.
- 2. Open the valve the liquid flows in.
- 3. Close the valve.
- 4. Pull out the dip sampler, take out the sample.



### Dip Sampler, plunging siphon

Volume ml	Immersion depth cm	Tube Ø mm	Item no.	
140	50	25	5315-0050 N4	
280	100	25	5315-0100 N4	
500	100 - 200	25	5315-0200 N4	
	Volume ml 140 280	140 50 280 100	Volume ml Immersion depth cm Tube Ø mm   140 50 25   280 100 25	Volume ml Immersion depth cm Tube Ø mm Item no.   140 50 25 5315-0050 N4   280 100 25 5315-0100 N4