Model MK 720 | Dynamic climate chambers for rapid temperature changes

The BINDER MK series chamber is suitable for all heat and cold tests between -40 °C and 180 °C. The MK is a clever alternative to complex individual solutions for cyclical temperature tests.

BENEFITS

- Homogeneous climate conditions thanks to APT.line™ technology
- · Comprehensive programming and data acquisition
- · Large heated viewing window



Model 720

MAIN FEATURES

- Temperature range: -40 °C to +180 °C
- APT.line™ preheating chamber technology
- Programmable condensation protection for test material
- · Heated viewing window with LED interior lighting
- BINDER Multi Management Software APT-COM™ Basic Edition
- Troubleshooting system with visual and audible alarms
- Intuitive touchscreen controller with time-segment and real-time programming
- Internal data logger, measured values can be read out in open format via
- 2 access ports with silicone plugs, 80 mm, left and right
- Class 2 independent adjustable temperature safety device (DIN 12880) with visual alarm
- · 4 castors, two with brakes
- Computer interface: Ethernet
- Adjustable ramp function
- Integrated chart recorder
- Real-time clock
- Door heating
- 1 stainless steel rack
- · Inner chamber made of stainless steel
- CFC-free refrigerant R-452A
- Cooling with compressor cooling unit



Model 720

OPTIONAL EQUIPMENT

- APT-COM Multi Management Software for managing, recording, and documenting unit parameters
- Racks option of stainless steel racks, depending on test requirements
- · Compressed air supply system
- Services wide range of services to ensure that the unit functions correctly

ORDERING INFORMATION


Interior volume [L]	Voltage	Option model	Version	ArtNo.
	400 V 3~ ph 50 Hz	Standard	MK720-400V	9020-0377
734	480 V 3~ ph 60 Hz	with voltage and frequency converter	MK720-400V MK720-480V-C	9020-0356

TECHNICAL DATA


Description	MK720-400V1	MK720-480V-C1
Article Number	9020-0377	9020-0356
Performance Data Temperature		
Temperature range [°C]	-40180	-40180
Temperature variation depending on setpoint [± K]	0.32	0.32
Temperature fluctuation depending on setpoint [± K]	0.10.5	0.10.5
Average heating-up rate according to IEC 60068-3-5 [K/min]	4	4
Average cooling down time according to IEC 60068-3-5 [K/min]	4.5	4.5
Max. heat compensation at 25 °C [W]	3500	3500
Electrical data		
Rated Voltage [V]	400	480
Power frequency [Hz]	50	60
Nominal power [kW]	8.7	8.7
Unit fuse [A]	16	16
Phase (Nominal voltage)	3~	3~
Measures		
Interior volume [L]	734	734
Net weight of the unit (empty) [kg]	570	570
Permitted load [kg]	160	160
Load per rack [kg]	40	40
Viewing window width [mm]	508	508
Viewing window height [mm]	300	300
Wall clearance back [mm]	300	300
Wall clearance sidewise [mm]	200	200
Internal Dimensions		
Interior width [mm]	1200	1200
Interior height [mm]	1020	1020
Interior depth [mm]	600	600
Doors		
Unit doors	1	1
Housing dimensions not incl. fittings and connections		
Width net [mm]	1580	1580
Height net [mm]	2005	2005
Depth net [mm]	1140	1140
Environment-specific data		
Energy consumption at 20 °C	1900 Wh/h	1900 Wh/h
Sound-pressure level [dB(A)]	65	67
Fixtures		
Number of shelves (std./max.)	1/11	1/11


¹ All technical data is specified for unloaded units with standard equipment at an ambient temperature of +22 °C ±3 °C and a power supply voltage fluctuation of ±10 %. The temperature data is determined in accordance to BINDER factory standard following DIN 12880, observing the recommended wall clearances of 10 % of the height, width, and depth of the inner chamber. Technical data refers to 100 % fan speed. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.

DIMENSIONS incl. fittings and connections [mm]

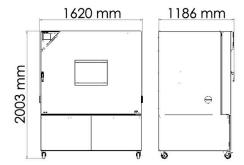

CHARTS

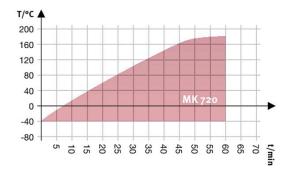

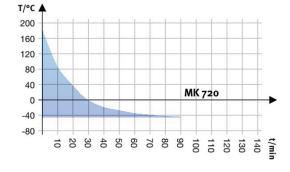

Heating up rate



Cooling down rate

Heat compensation chart

OPTIONS

Designation	Description	*	ArtNo.
	left		
	30 mm	01	8012-1322
	50 mm	01	8012-1328
	8o mm	01	8012-1334
	100 mm	01	8012-1542
	125 mm	01	8012-1351
	right		
Access port with silicone	30 mm	01	8012-1319
olug	50 mm	01	8012-1325
	8o mm	01	8012-1331
	100 mm	01	8012-1539
	125 mm	01	8012-1348
	top		
	8o mm	01	8012-1536
	100 mm	01	8012-1530
	125 mm	01	8012-1533
Analog output 4-20 mA	for temperature values (output not adjustable)	-	8012-108
Calibration certificate, expanded	for temperature; for extending the measurement in center of chamber to include another test temperature	-	8012-1124
	for temperature, measurement in center of chamber at specified temperature	_	8012-1143
Calibration certificate,	temperature measurement incl. certificate and 27 measuring points at specified temperature	-	8012-160
emperature	temperature measurement incl. certificate, 15- 18 measuring points at specified temperature	-	8012-1581
	temperature measurement incl. certificate, 9 measuring points at specified temperature	-	8012-1560
Compressed air	for the connection to an existing pressurized air network	-	8012-1797
Door lock	lockable door handle	-	8012-186
Ory-air purge	uncontrolled, incl. connection; for preventing condensation	-	8012-186
Notch-type access port	notch-type access port in door, 100 x 35 mm	-	8012-1851
Pt 100 temperature sensor	additional flexible Pt 100, interior, for displaying the temperature on the unit display	-	8012-109
RS 485 interface, 2-wire	Additional serial interface can be used parallel to Ethernet, for Multi Management Software APT-COM™	-	8012-1768
Shelf, perforated	Stainless steel	-	8012-225
Temperature safety device class 2	with visual alarm (DIN 12880)	-	8012-1792
Zero voltage relay contacts	for controlling 3 relay contacts via program regulators, accessible via 6-pin DIN socket (max. 24 V - 2.5 A)	-	8012-109
Notes > See last page			

^{*} Notes > See last page

ACCESSORIES

Designation	Description	*	ArtNo.
APT-COM™ 4 GLP- Edition	for working under GLP-compliant conditions. Measured values are documented in a tamper-proof way in line with the requirements of FDA Regulation 21 CFR 11.		
	version 4, GLP edition	-	9053-0042
APT-COM™ 4 PROFESSIONAL-Edition	convenient unit and user management built on the BASIC edition. Suitable for networking up to 100 units.		
	version 4, PROFESSIONAL edition	-	9053-0040
pH-neutral detergent	concentrated, for gentle remove of residual contaminants; 1 kg	-	8012-2250

^{*} Notes > See last page

Designation	Description	*	ArtNo.
	IQ/OQ documents – supporting documents for validation performed by customers, consisting of: IQ/OQ checklists incl. calibration guide and comprehensive unit documentation; parameters: temperature, CO ₂ , O ₂ , pressure, depending on unit		
	Digital in PDF format	-	7057-0001
Dualification de sumante	Hard copy inside folder	-	7007-0001
Qualification documents	$IQ/OQ/PQ$ documents – supporting documents for validation performed by customers, according to customer requirements, PQ section added to qualification folder IQ/OQ ; parameters: temperature, CO_2 , O_2 – or pressure, depending on unit		
	Digital in PDF format	-	7057-0005
	Hard copy inside folder	-	7007-0005
Rack	stainless steel	-	8012-2054
Rack accessories	fasteners (1 set of 4) for additional security of racks	-	8012-2280
Rack, reinforced	stainless steel, with fasteners (1 set of 4)	-	8012-0684
	RS 422 cable set and RS 485 / RS 422 interface converter for connection to 10-way plug distributor		
RS 485 / RS 422 nterface converter	115 V option model	-	8012-0599
interiace converter	230 V option model	-	8012-0589
Shelf, perforated	Stainless steel	-	8012-2253
The kit contains a client bridge that establishes a wireless connection between BINDER units and APT-COM4, LIMS, or customer-specific software via an Ethernet interface. It provides an alternative solution in situations where units need to be positioned in locations without an on-site Ethernet connection. In secured networks, installation and configuration must be performed by the customer's IT service.		-	8012-2262

^{*} Notes > See last page

SERVICES

Designation	Description	*	ArtNo.
Maintenance contracts			
Calibration services			
Validation services			
Installation services			
Maintenance services			
Warranty service			
1-year warranty extension	The warranty is extended by 1 year from the delivery date, wear parts are excluded	-	DL50-0040
BRONZE 3-year maintenance contract	Maintenance service as contractually agreed, visual inspection of mechanical and electrical components, check of control response, 20% discount on spare parts	05	DL20-0710
Calibratian and Grant American	Calibration of one (1) test temperature specified by the user in center of chamber, including certificate	03, 04, 05	DL30-0101
Calibration certificate, temperature	Extension of calibration of one (1) additional test temperature specified by the user in the center of the usable space, including certificate	03, 04, 05	DL30-0102
Execution of IQ/OQ	Execution of IQ/OQ in accordance with qualification folder	05	DL41-0200
Execution of IQ/OQ/PQ	Execution of IQ/OQ/PQ in accordance with qualification folder	05	DL44-0500
GOLD 3-year maintenance contract	Maintenance service as contractually agreed, visual inspection of mechanical and electrical components, check of control response, 20% discount on spare parts, testing of all key functions, replacement of wear parts, calibration of one temperature and CO_2 value, including certificate	05	DL20-0940
Leakage test	Testing and logging of the leak-tightness of systems containing hydrofluorocarbons in accordance with Regulation EC842/2006, incl. test book	05	DL00-0034
Maintenance	One-off maintenance service in accordance with maintenance schedule. Visual inspection of mechanical and electrical components, testing of all key functions. Calibration of a test temperature specified by the user in center of usable space without certificate	05	DL20-0500
SILVER 3-year maintenance contract	Maintenance service as contractually agreed, visual inspection of mechanical and electrical components, check of control response, 20% discount on spare parts, testing of all key functions, calibration of one test temperature specified by the user in the center of the usable space, without certificate	05	DL20-0830
Temperature measurement, 18 measuring points	Temperature measurement with 18 measuring points with a set value specified by the user, including certificate	03, 04, 05	DL30-0118

^{*} Notes > See last page

Data Sheet Model MK 720

Designation	Description	*	ArtNo.
Temperature measurement, 27 measuring points	Temperature measurement with 27 measuring points with a set value specified by the user, including certificate	03, 04, 05	DL30-0127
Temperature measurement, 9 measuring points	Temperature measurement with 9 measuring points with a set value specified by the user, including certificate	03, 04, 05	DL30-0109
Connect the unit to the customer-side connections (electricity, water, wastewater, gas), basic functions Unit commissioning check, brief operating instructions. (excl.: unpacking, setup, controller instructions, programming, installation work)		05	DL10-0300
Unit instructions	Instruction regarding operating principle and basic functions of the unit, operation of the control electronics including programming	05	DL10-0700

^{*} Notes > See last page

NOTES

- Condensation may occur in the area around the access port. Access ports may be placed in custom locations for an additional charge. UL mark is not granted when this option is used.
- 02
- Sensor calibration is performed in an accredited calibration laboratory. 03
- Calibration is performed according to the BINDER factory standard.
- Quoted prices do not include travel costs. Please refer to the chapter on BINDER Service for travel costs for your region. Quoted prices for services performed in Switzerland do not include a country-specific added fee (available on request).

BINDER GmbH

Tuttlingen, Germany TEL +49 7462 2005 0 FAX +49 7462 2005 100 info@binder-world.com www.binder-world.com

BINDER Asia Pacific (Hong Kong) Ltd.

Kowloon, Hong Kong, P.R. China TEL +852 39070500 FAX +852 39070507 asia@binder-world.com www.binder-world.com

BINDER Environmental Testing Equipment (Shanghai) Co., Ltd. Shanghai, P.R. China TEL +86 21 685 808 25 FAX +86 21 685 808 29 china@binder-world.com www.binder-world.com

BINDER Inc. Bohemia, NY, USA TEL +1 631 224 4340 FAX +1 631 224 4354 usa@binder-world.com www.binder-world.us